Pipette Filter Plugs Art. No. 100220

Filter Plugs protect the pipette and the sample from contamination by any fluids, for example by accidental splashes or droplets during pipetting, foaming samples, etc. They reduce the service interval and increase lifetime of the pipette. The strongly hydrophilic, porous filter material is a highly efficient absorber and acts as an efficient barrier to prevent sample components from attaching to the pipette cone or entering into the pipette. Filter Plugs are the cost-effective alternative to filter tips for many applications. The universal Filter Plugs fit into 1 ml air displacement pipettes with a tip cone opening of 3 mm (-) 4.5 mm like: Eppendorf, Biohit/ Sortorius, Brand, Capp, Thermo Finn pipette, etc.


Roll the Filter Plug lengthwise between thumb and forefinger.

Note: Do not roll too powerful, so that the filter structure is not damaged. Disinfect hands or wear gloves.

Completely insert the rolled Filter Plug into the pipette. The Filter Plug now expands in the pipette. Wait for 30 – 60 seconds until the Filter Plug is stucked.

Recommendation: Before inserting a Filter Plug clean tip cone opening for example with 70% ethanol.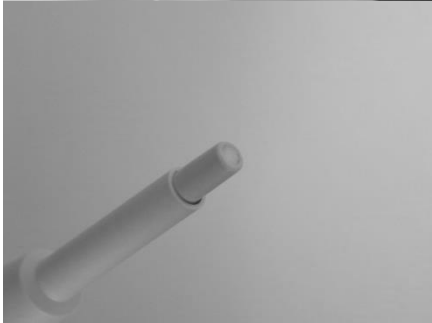

Insert Filter Plug as flush as possible with the tip cone opening. For some pipettes the Filter Plug may be too long. If necessary, cut off the protruding end or shorten the Filter Plug prior to insertion.

Note: Before usage clean scissors, scalpel, etc. for example with 70% ethanol.

For exchange pull the used Filter Plug with a needle, cannula, pointed tweezers, etc. out of the pipette.

The interval of changing the Filter Plugs depends on the application and the samples: after about 500 pipettings, visible contaminations of the white filter surface or determined over-aspiration.

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